PHENOLPHTHALEIN TEST KİTİ

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PHENOLPHTHALEIN TEST KIT

Phenolphthalein is a chemistry widely used in laboratory and field studies to show the presence of blood. Phenolftalein reacts with the heme group in blood and gives pink color.

With this test kit, blood stains can be determined easily and quickly using the phenolphthalein method.

CONTENT OF TEST KIT	
Alcohol Reagent	30 mL
Phenolphthalein Reagent	30 mL
Hydrogen Peroxide Reagent	30 mL
Distilled Water –15 mL	
Swab – 50 Pcs	
Filter Paper – 50 Pcs	
Example of known Blood Stain - 1 Pc	

PHENOLPHTALEIN TEST METHOD

❏ Sampling:

A drop of **Alcohol Reagent** is placed on a cotton swab or filter paper and then rubbed onto a stained surface.

❏ PROCEDURE:

Add 1 drop of **Phenolphthalein Reagent**, and 1 drop of **Hydrogen Peroxide Reagent** onto the sampled cotton swab or filter paper.

❏ RESULT:

- **Positive Test:** If the color of the cotton swab or filter paper turns pink immediately, the result is interpreted as **POSITIVE TEST**. So the sample is blood.
- **Negative Test:** If the color of the cotton swab or filter paper lasts for 1 minute or longer, the result is interpreted as **NEGATIVE TEST**. So the sample is not blood.
- You can observe a positive result by experimenting with a known blood stain sample before performing the experiment.